## Goniometer Goniometers

MOGO 65-40-65

/ 23.11.2020

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**Users are required to follow the safety and security advice.**

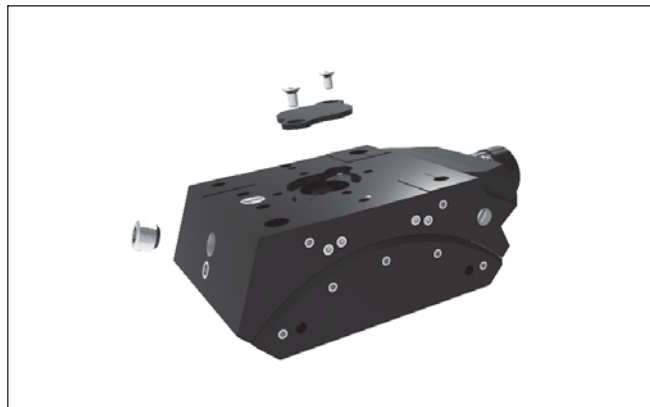
Lubricating greases consist primarily of a base oil and a thickener. Not every base oil and not every thickener can be mixed with each other. When combining immiscible components, you get an inhomogeneous mixture that may have a low dropping point or may tend to harden. Thus, mixtures of different lubricating greases should be avoided. It is highly recommended that the original lubricating greases are used for relubricating.

The **worm gear** and the **guides** are lubricated by us with the **OWIS® SST.F8** grease.

Lubricating grease SST.F8 in a 5 ml applicator, part no. 90.999.0008

Relubrication should take place after about 250 operation hours, at least every 3 months.

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1. Turn off the control unit and make sure that it cannot be turned on accidentally.
2. Remove both screws of the bottom cover plate and the screw of the lateral bore.  
**Attention!** During the actual lubrication procedure the MOGO must be halted, so that the applicator needle cannot be seized and pulled in by the worm.
3. Apply a thin thin film of lubricating grease SST.F8 through the bottom opening.
4. Insert the applicator needle into the open bore up to the worm and fill in about 0.1 ml grease SST.F8 (corresponds to ½ applicator scale-line).
5. Turn the gonio another quarter of the travel and fill again about 0.1 ml grease SST.F8 into the opening. Continue this procedure about the complete travel range of the gonio.
6. Apply evenly a thin film of SST.F8 on the guides.
7. Fix the screw of the lateral bore and the bottom cover plate.