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Using adjustable inserts with fine-thread screws JT 65F and with disc springs JT 65T in conjunction with holders for mounted optics OH 65-D45 or cubes W 65, highly stable beam deflectors can be assembled. With cubes W 65, it is a simple matter to build a microscope setup with two viewing points. By combining several cubes, precise ring resonators, Mach-Zehnder, Michelson interferometers, etc. can be built.

The adjustable inserts with fine-thread screws JT 65F can be set without the need of any tools. They are especially suitable for setups where frequent adjustment is necessary. Two fine-thread screws permit optical elements to be tilted in two axes. A third screw enables an additional displacement of the centre of rotation.

In the adjustable inserts with disc springs JT 65T the optic holders are supported on packets of 3 disc springs against the flange component, and can be adjusted. They are very precise, shock and vibration resistant and long-term stable. Also here two 2 packets enables the adjustment of the optics in theta and phi. The third screw provide an additional displacement of the centre of rotation.

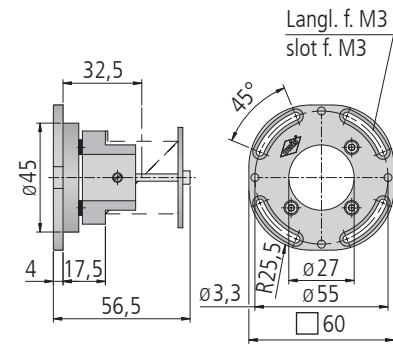
These adjustments allow compensation of different thickness, or the setting of beamsplitter cubes at a defined inclination of the front and exit field, to eliminate reflections.

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Adjustable Inserts for Beamsplitter Cubes and Prisms

with 30 mm side. By the use of lateral clamping, 30 mm prisms can also be installed.

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
Adjustable Inserts for Beamsplitter Plates

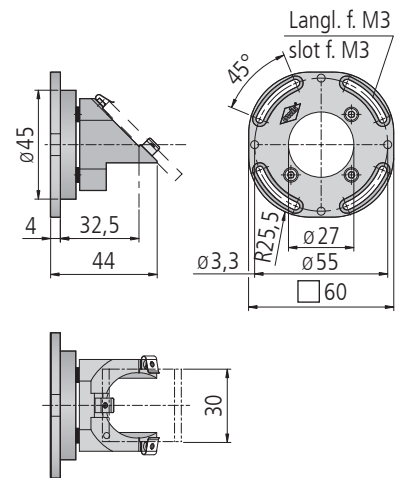with 30 x 42.5 x 4 mm. The plates are retained by springs at three points. The low-stress mounting means that the insert can be used for interferometry, even at large beam diameters.

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
Adjustable Inserts for Deflection Mirrors

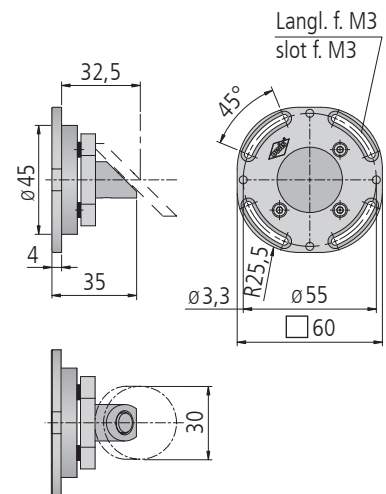with 30 x 42.5 x 4 mm for deflecting beams at 90°. Low-stress bonding to the optical glass substrate makes this suitable for high-quality laser mirrors. Thermal stresses are absorbed in the substrate, and so do not affect the mirror. The optimum mirror thickness is 4 mm.

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
Adjustable Inserts for Front Surface plane and concave Mirrors

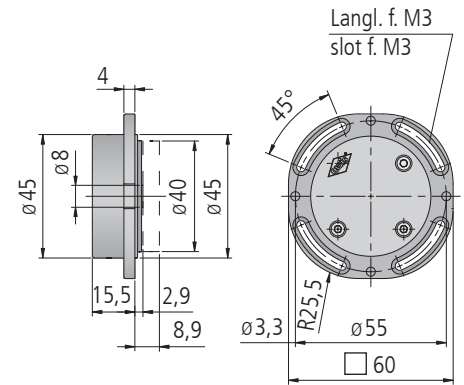with diameter 40 mm and plane backside. Low-stress bonding of the mirror to the optical glass substrate makes this suitable for interferometry.

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
Spacer Ring 9.5 mm

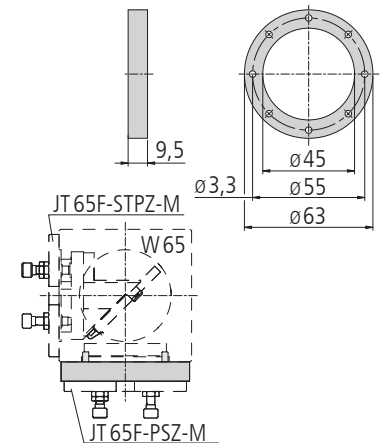when fitting two adjustable inserts to one cube W65, for example a beamsplitter and a plane mirror, the 9.5 mm spacer ring is required.

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