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Users are required to follow the safety and security advice.

Lubricating greases consist primarily of a base oil and a thickener. Not every base oil and not every thickener can be mixed with each other. When combining immiscible components, you get an inhomogeneous mixture that may have a low dropping point or may tend to harden. Thus, mixtures of different lubricating greases should be avoided. It is highly recommended that the original lubricating greases are used for relubricating.

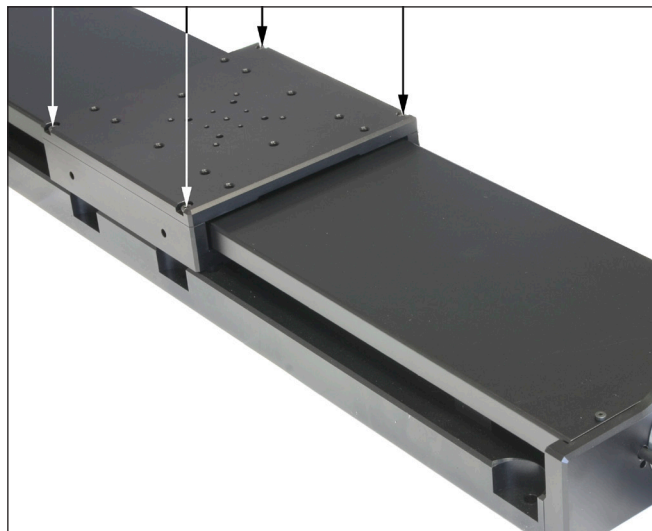
The **spindle** resp. the **ball screw** are lubricated by us with the **OWIS® SST.F11** grease.
Lubricating grease SST.F11 in a 5 ml applicator, part no. 90.999.0011

The **recirculating ball bearing guides** are lubricated by us with the **OWIS® SST.F2** grease.
Lubricating grease SST.F2 in a 5 ml applicator, part no. 90.999.0002

Relubrication should take place after about 500 operation hours or after 100 km total travel, at least every 3 months.

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
Lubrication procedure


1. Turn off the control unit and make sure that it cannot be turned on accidentally.
2. Loosen the fixing screws of the mounting plate, after (if necessary) removing the setup. The mounting plate is pinned. Therefore, please take it off perpendicularly, so that the fitting accuracy is not affected.

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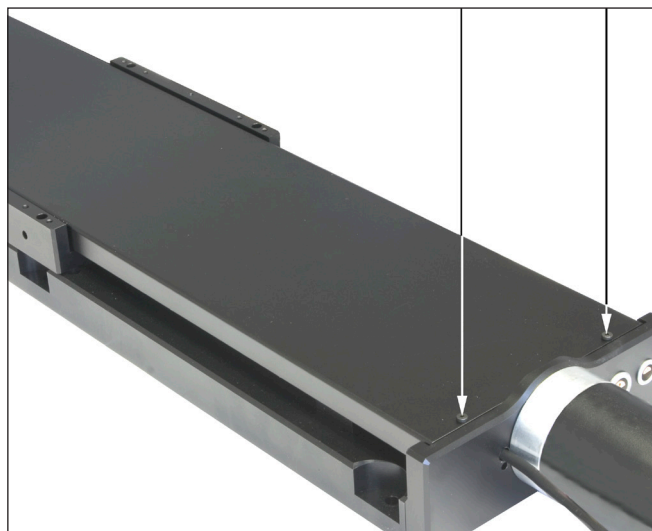3. Screw off the metal covering and slide it out backward.

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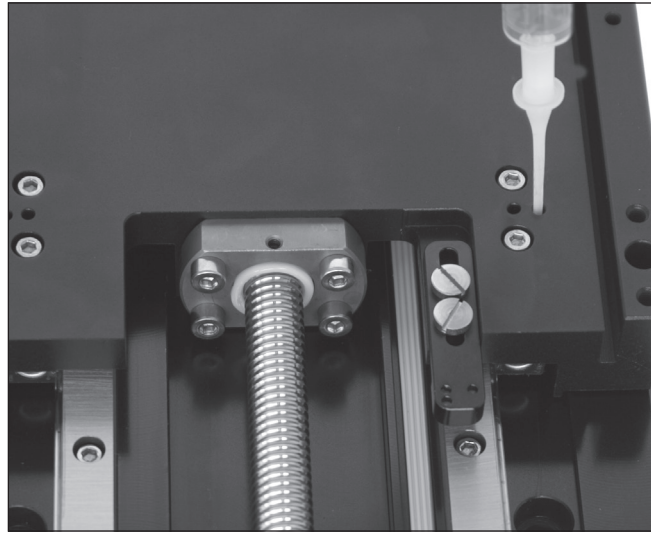4. Remove the used grease or dust particles from the guides and the threads of the spindle resp. the ball screw with a dry, lint-free cloth. Cleansers such as benzine or petroleum are normally not necessary.
5. Each of the four recirculating ball bearing slides has a hole where the lubricating grease SST.F2 has to be injected by means of an appropriate syringe.

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6. Apply evenly a thin film of lubricating grease SST.F11 onto the thread surface. Dispense the grease with a brush carefully on the complete surface.

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7. Assemble the linear stage in reverse order.